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Dialog box titled "Edit Cell Type". It contains fields for Cell Line, Name, Cell Type 1, Cell Type, Lot, Source, Prepared By, Prepared On, Passage, and Thawed. The "Seeding Density" field is highlighted with a red box.

The image shows two overlapping windows. The top window is the "Protocol" screen, with the "Custom" tab selected. It shows "3 Measurement Cycles" and "0 Commands". The "0 Commands" section is highlighted with a red box. The bottom window is the "Assay Summary" screen, showing "Project Information", "Plate Information", and "Notes". A red arrow points from the "0 Commands" section to a warning message on the right side of the Assay Summary screen.

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Plate Map

The Plate Map interface shows a 12x8 grid of wells. Wells D5 and E8 are highlighted with a red box. The interface includes buttons for "Clear Plate" and "Distribute Groups".

Note: Well D5 and E8 must be seeded with cells for use with Cytation and the Cell Counting and Normalization Application.
Select a group from the left, then click or drag with the mouse to apply that group to the plate map.

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